Our knowledge - Your success

Bedienungsanleitung / Instructions / Mode d'emploi / Manual

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The control unit HT 50 is used in combination with a heated microscope stage, a heated laboratory plate or a heated storage container.

Equipment:

The control unit HT 50 is equipped with an electronic two-step-controller with digital LED display. The controller is operated by a membrane keyboard.

Instructions

Control unit HT 50

Setting into operation:

The device should be used in dry rooms only.

The control unit must be adjusted to the corresponding voltage (230 V or 115 V). The actual voltage setting may be read and/or adjusted at the integrated fuse holder of the power cord connector at the rear side. In order to change these settings it is necessary to pull out the fuse holder from the casing. Then pull out the fuse socket from the fuse holder. Rotate end to end and replace it so that the proper voltage is displayed through the window. Please do not forget to insert the proper fuse, before putting back the fuse holder into the casing. (Please watch the value of the fuses when changing the input voltage!!)

IMPORTANT: Use only fuses **315 mA(T)** DIN 41661 with 230 V / 50 Hz,
and only fuses **630 mA(T)** DIN 41661 with 115 V / 60 Hz

Connect the respective plug of the external heating system to the unit and switch on. The output jack is performance-oriented assigned to the heating system. Do never mix up control units and heating systems. The „ON/OFF“ switch is at the rear side of the unit. After switching on, the controller starts automatically a self-test, then the actual value is indicated.

Adjustment of the set value

When pushing the „SET“ button once quickly, the set value is indicated for about 5 seconds.

Change of the set value:

Press the „SET“ button for at least 2 seconds. After that, the set value is indicated, while a LED above the decimal point flashes. The actual temperature value may now be adjusted with the buttons „▲“ or „▼“. The desired set value is automatically stored after 15 seconds without pushing any button respective after having pushed once the „SET“ button. After that, the actual value is indicated again.

Your contact for further technical inquiries:

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GmbH & Co. KG
EMV-Testhaus GmbH
Minitüb Abfüll- und Labortechnik Verwaltungs GmbH